



T CELL RESTIMULATION

OVERVIEW

Cell cultures showing signs of exhaustion (typically at day 7–10 of expansion) can be restimulated several times by adding fresh CD3/CD28 Macrobeads™ (cat. # PM-CAR2002, ProMab Biotechnologies) and recombinant IL-2 (cat. # Pr21269, ProMab Biotechnologies). The CD8+ T cells remain cytotoxic after repeated restimulations. Restimulation is typically necessary when cell shrinking, and a reduced rate of proliferation are observed.

MATERIALS NEEDED

1. Buffer: Phosphate buffered saline with 0.1% bovine serum albumin and 2 mM EDTA, pH 7.4 (PBS w/0.1% BSA).
2. CD3/CD28 Macrobeads™ (cat. # PM-CAR2002, ProMab Biotechnologies).
3. Magnet
4. Culture medium: Advanced RPMI Medium 1640 with 2 mM L-Glutamine, 10% FCS/FBS and 100 U/ml penicillin/streptomycin can be used. Alternatively, Cancer Stem Premium™ (cat. # 20101, ProMab Biotechnologies) with 100 U/ml penicillin/streptomycin, or another equivalent culture medium.
5. Recombinant human IL-2 (cat. # Pr21269, ProMab Biotechnologies).
6. Heat inactivated Fetal Calf Serum (FCS).
7. Flat bottom tissue culture plates or tissue culture flasks.
8. Humidified CO₂ incubator.

WASHING OF MACROBEADS™ BEFORE USE

1. Resuspend the Macrobeads™ in the vial (i.e. vortex for >30 sec, or tilt and rotate for 5 min).
2. Transfer the desired volume of Macrobeads™ to a tube.
3. Add an equal volume of buffer, or at least 1 mL, and mix (vortex for 5 sec, or keep on a roller for at least 5 min).
4. Place the tube on a magnet for 1 min and discard the supernatant.
5. Remove the tube from the magnet and resuspend the washed Macrobeads™ in the same volume of culture medium as the initial volume of Macrobeads™ taken from the vial.

Additional Products and Services:

- [Mouse Monoclonal Antibody](#)
- [Rat Monoclonal Antibody](#)
- [Rabbit Monoclonal Antibody](#)
- [Human Monoclonal Antibody](#)
- [Polyclonal Antibody](#)
- [Antibody Sequencing](#)
- [Hybridoma Sequencing](#)
- [CAR T-cells](#)
- [Lentivirus production](#)
- [Cancer Stem Cells](#)
- [Specialty Cell Culture Media](#)
- [T-cell Expansion beads](#)

Ask about our full line of CRO services to provide supplemental assistance or the entire support necessary to complete your project on time and with the data you need to move forward.



T CELL RESTIMULATION



RESTIMULATION

1. Prior to restimulation, remove the used Macrobeads™ by transferring the cells to a suitable tube.
 2. Place the tube in the magnet for 1–2 min.
 3. Transfer the supernatant containing the cells to a new tube.
 4. Split the cultures back to a density of $0.5\text{--}1 \times 10^6$ cells/mL in culture medium containing 300 U/mL rIL-2 and repeat the Expansion procedure.
- Guidelines for restimulation are provided in Table 1. Optimize for your application. Do not use an excess volume of Macrobeads™, as excess Macrobeads™ may inhibit expansion.

RESTIMULATION GUIDELINES

Number of Cells	1×10^4 T-Cells
Cell type	Subsequent restimulations*
CD4 ⁺ (polyclonal)	8-10 day intervals
CD8 ⁺ (polyclonal)	7-9 day intervals
T cells	7-9 day intervals*

Table 1. Restimulation guidelines for anti-CD3/CD28-expanded cultures



T CELL ACTIVATION AND EXPANSION



BEAD-TO-CELL RATIO

Type of culture plate/flask	24-Well Plate	175 cm ² Tissue Culture Flask
Cell concentration	1 × 10 ⁶ T cells/well	50 × 10 ⁶ T cells/flask
MacroBeads™	25 μL	1,250 μL
rIL-2	300 U/mL	300 U/mL
Seeding volume (medium)	1–2 mL	50–100 mL

Table 2. Volume recommendations for bead-to-cell ratio = 1:1